

Shortcuts for Vicon Nexus

- F1 ----- Vicon Nexus help
- F2 ----- Data Management
- F4 ----- Quick Report window
- F5 ----- Full screen
- F6 ----- Sounds dialog box
- F7 ----- Options dialog box
- F8 ----- System Preparation Tools pane
- F9 ----- Subject Preparation Tools pane
- F10 ----- Capture Tools pane
- F11 ----- Label/Edit Tools pane
- F12 ----- Pipeline Tools pane


Useful hotkeys

- Ctrl+Enter ----- Start/Stop capture
- Ctrl+Tab ----- Switch Live/Offline mode
- Ctrl+Space ----- Display marker labels
- Ctrl+← ----- Move to previous event
- Ctrl+→ ----- Move to next event
- Ctrl+Z ----- Undo
- Ctrl+Y ----- Redo
- Ctrl+S ----- Save trial
- Ctrl+R ----- Reset Core Processor
- Space ----- Pause/Restart real-time data streaming
- Space ----- Play/Stop offline data
- Esc ----- Exit current mode (labeling, etc.)
- Ctrl+Shift+B ----- Reset bumped status of all cams

Mouse Actions

- | | |
|-------------------------------------|--------------------------------------|
| Right mouse click and drag | Zoom 3D-Space |
| Left mouse click and drag | Rotate 3D-Space |
| Left and right mouse click and drag | Move3D-Space |
| Left mouse click | Select individual item |
| CTRL+Left mouse click | Select several items |
| ALT + Left mouse click and drag | Select items inside the bounding box |

CGM Marker Placement

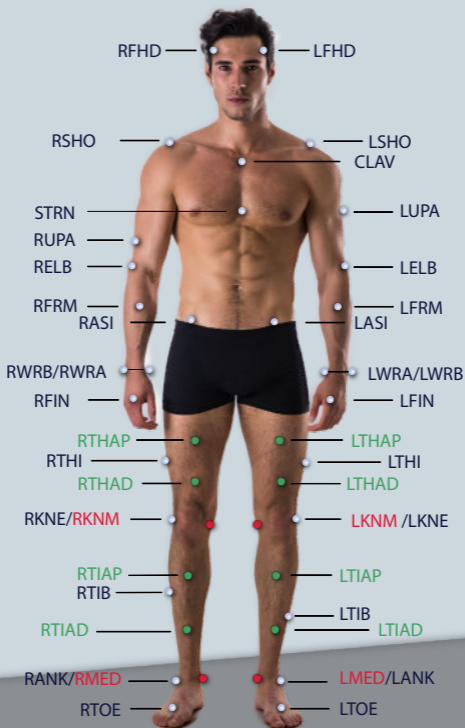

● PiG ● CGM 2.1 + ● CGM 2.3

